


## Hei-VOLUME Distimatic 24/7 Industrial with adaptor for A-series glassware

P/N: 036064600

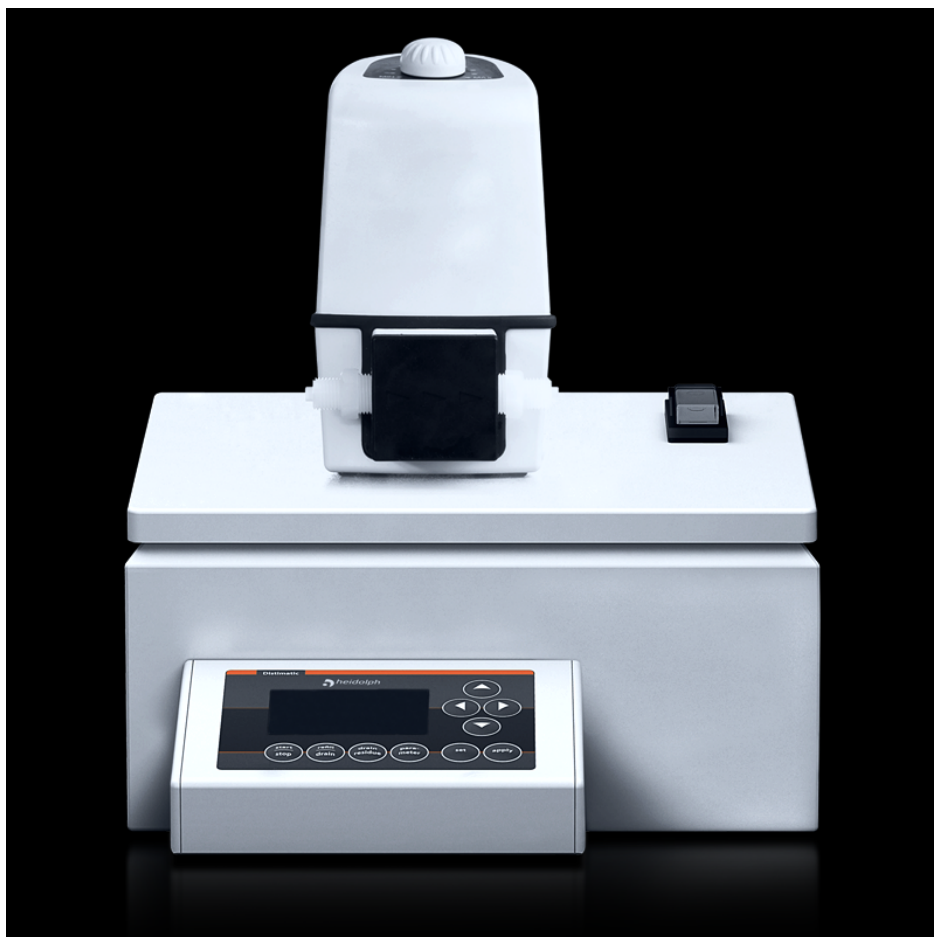

For high-performance distillation with the best separation performance, suitable for most solvents (A1 version), also when processing constantly foaming media (A version) or for foam suppression at the start of the process (A2 version): Version with adapter for Hei-VAP Industrial with glassware of the A series (A, A1, A2)

You can use the control panel on the control box of the Hei-VOLUME Distimatic 24/7 module to enter the parameters for the sensor or time-controlled filling of the rotary flask. The emptying times of the coated collector vessel (approx. 700 ml) and the residue from the rotary flask by the two chemical membrane pumps can also be programmed. The control box can be mounted directly on a Hei-VAP Industrial large-scale rotary evaporator or on a nearby wall using the supplied bracket with hose guide.

In combination with the corresponding evaporator system, the Hei-VOLUME Distimatic 24/7 automatic module enables efficient, unattended operation around the clock for every range of application.

For the greatest possible safety and a wide range of applications, all parts that come into contact with the media, including the supplied tubing, are made of chemical-resistant materials.



## Hei-VOLUME Distimatic 24/7 Industrial - Technical Data

|                                   |   |
|-----------------------------------|---|
| Permissible ambient conditions    | 5 – 31 °C at 80 % rel. humidity 32 – 40 °C decreasing lineary up to max 50% rel. humidity |
| Weight                            | 5 kg  |
| Protection class (EN 60529)       | IP 65 (control panel and residue pump) IP 54 (Housing)                                    |
| Display                           | Separate control panel; digital   |
| Power input                       | 3500  |
| Dimensions (w/d/h)                | 300 x 200 x 120 mm (approx.; without mounted components)                                  |
| Compatibility                     | Hei-VAP Industrial and Laborota 20 series   |
| Automatic refill of sample        | yes   |
| Automatic discharge of condensate | yes   |
| Automatic drainage of residue     | yes   |
| Emergency-Stop button             | yes   |
| Functional Principle              | internal (condensate pump) and external diaphragm pump (residue pump)                     |